Manual Cryo Labels | Kryo-Etiketten

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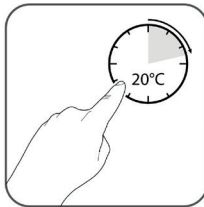
Right handling of the Cryo Labels


Temperature Resistance

The cryogenic labels withstand temperatures from minus 196 °C to plus 121 °C and are suitable for long-term storage in liquid nitrogen and transport on dry ice. Even after repeated freezing and thawing, the cryogenic labels have shown good adhesion. If a label partly peeling off, it is possible to fix it again by pressing. The cryo-labels can be attached to a frozen tube. You should nevertheless dry the surface thoroughly first.

Applying the Labels

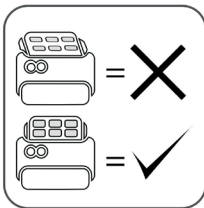Make sure that the surface of your vessel or tube is not damp before applying the label. To prevent the glue from drying out, apply the label at room temperature immediately after removing it from the backing film. For getting the best result, labels should be fixed at least 30 minutes before application at room temperature (approx. 20°C). If possible, stick the labels slightly overlapping.

Print Information

Please make sure that the label sheets do not slip when printing, to avoid any damage. Use the enclosed adjustment template to align your printer.

Chemical Resistance

Materials and Methods

Cryo Labels, format 28×12 mm/ white, marked with:

- Laser printing (HP Color Laser Jet CM 1312),
- Ball-point pen (no-name product),
- Pencil (Rheita, HB)

Labels were placed on 2 ml tubes and submersed in the respective solvent for 24 h. Shorter exposition times may increase resistance.

Overview Chemical Resistance at Room Temperature

Solvents	Adhesion	Laser printing		Ball-point pen		Pencil	
		Legibility	abrasion resistance	Legibility	abrasion resistance	Legibility	abrasion resistance
Acetic acid 25 %	G	G	G	G	G	G	L
Acetone	N	L	L	L	L	G	L
Acetonitrile	L	G	G	L	G	G	L
Ammoniumhydroxide (1M)	G	G	G	G	G	G	L
Ammoniumhydroxide (25%)	G	G	G	L	G	G	L
Ethanol	G	G	G	G	G	G	L
Ethanol 70 %	G	G	G	G	G	G	L
Formic acid (25%)	G	G	G	L	G	G	L
Hexane	N	G	G	G	G	G	L
Hydrochloric acid (10%)	G	G	G	G	G	G	L
Hydrogenperoxide (30%)	G	G	G	G	G	G	L
1-Propanol	L	G	G	G	G	G	L
Methanol	G	G	G	G	G	G	L
Methylenchloride	N	L	L	N	N	G	L
Phosphoric acid (25%)	G	G	G	G	G	G	L
Potassium Hydroxide (KOH, 1M)	L	G	G	G	G	G	L
Potassium Hydroxide (KOH, 32%)	L	L	L	L	L	L	L
Sodium hydroxide (1M)	G	G	G	G	G	G	L
Tetrahydrofurane (THF)	N	N	N	G	G	G	L
Toluene	N	L	N	G	G	G	L
Water distilled deionized	G	G	G	G	G	G	L

G	Good chemical resistance
L	Limited chemical resistance
N	No chemical resistance, use not recommended

Use of the Calibration Template with Microsoft Word template

You can download the Microsoft Word templates for the labels on the product pages of the scienova webshop.

- First open the template document in Microsoft Word.
- Enter a sample text and/ or illustration in one of the fields. Place the elements on the labels at least 1-1.5 mm from the edge.
- Advice: You can show the grid lines in the table properties . Click on the Table Layout tab and then under Settings on "show Gridlines". This will simplify the orientation and the optimal alignment of your printer to the template.
- Insert the supplied calibration template into the printer and print the labels. Attention: The positioning aids in the corresponding compartment must be set as precisely as possible to A4 format.
- If the printed frame differs from the frame line of the template, measure the horizontal and vertical displacement. This allows you to set a small offset in the settings menu of your printer (directly on the device).
- Repeat the process with the next table field until you have the best accuracy you can achieve.
- In Microsoft Word, you can move the template to the side and align it even more precisely: In the table tools under "Layout" → "Table Properties", you can easily move the entire table on the sheet. Test the different alignment variants [left, centered and right] and approach the optimal alignment.

Microsoft Windows, and Microsoft Word are trademarks of the Microsoft group of companies.

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scienova GmbH